Ion exchangers for demineralising water. Safe, reliable, top quality that can be regenerated time and time again.





# Ion exchangers from stakpure. The economic answer to demineralised water.



Whether for ultrapure water systems, automatic analysers, dishwashers, autoclaves – demineralised water is needed on a daily basis for all sorts of different purposes in medical practices, hospitals and laboratories. Whether in small quantities or up to 3000 litres an hour – we have the right ion exchanger for any need.

Inside a container, water flows over a mixed-bed of special resins that removes its salts and other minerals. A tried and proven method that guarantees water of a constant high quality all the time. The resins come with a long useful life and can be regenerated whenever necessary without any impact on the environment.

# For small quantities: Ion exchangers in cartridges of synthetic material.

These ion exchangers have been designed specifically for medical practices and small laboratories requiring up to 10 litres a day. The non-pressurised cartridge will be simply regenerated or replaced after a prolonged period of service. It is equipped with conductivity meter, set of hoses and wall mounting is optionally available.

### **Operating safety**

An innovative filter system prevents any resin from escaping. High product safety for downstream equipment as well.

stakpure ion exchangers	DS 450*	DS 750*
Model	one-way cartridge	regenerably cartridge
Flow rate	50 l/h	100 l/h
Capacity for 10° total dissolved solids	425 l	750 l
Pure water quality	0.1–20 µS/cm	0.1–20 µS/cm
Max. water temperature	30°C	30°C
Power supply	230 V/50-60 Hz	230 V/50-60 Hz
Material	Polyethylene	Polyethylene
Dimensions mm (Ø x height)	100 x 600	175 x 470
Connection	R ¾"	R 3⁄4″
Weight, approx.	3 kg	6 kg
Order no.	11042500	11170500
Replacement cartridge	11042700**	11175000
Wall mount	– included	13140400



\* conductivity meter included

\*\* set 2 pcs.

# For large quantities: lon exchangers in stainless steel cartridges.

For any situation requiring 300 to 3000\* litres of demineralised water an hour. A sophisticated flow system for incoming water ensures a fast and effective flow of water through the mixed-bed system with ion-exchange resins.

\* We recommend our reverse-osmosis systems in applications requiring higher quantities or where particular demands are placed on pure water.


### stakpure DS ion exchange cartridges

For up to 300 and 1000 l an hour. Stainless-steel cartridge. Hard-rubber carrying handles.



DS 1500	DS 2000	DS 2800	DS 4000	DS 6000
pressure-resistant				
pressure-resistant to 10 bar				
300 l/h	300 l/h	950 l/h	1000 l/h	1000 l/h
1500 l	2000 l	2800 l	4000 l	6000 l
0.1–20 µS/cm				
30°C				
Stainless steel – AISI 316				
240 x 410	240 x 490	240 x 570	240 x 700	240 x 1155
		R ¾ ″		
14 kg	18 kg	24 kg	27 kg	45 kg
12115000	12200000	12280000	12400000	12600000
13140500				
	300 l/h 1500 l 0.1–20 μS/cm 240 x 410 14 kg	pr 300 l/h 300 l/h 1500 l 2000 l 0.1–20 μS/cm 240 x 410 240 x 490 14 kg 18 kg	pressure-resistant   pressure-resistant to 10   300 l/h 300 l/h   300 l/h 300 l/h   1500 l 2000 l   0.1 – 20 µS/cm 30 °C   Stainless steel – AISI 31   240 x 410 240 x 490   240 x 410 240 x 490   14 kg 18 kg   12115000 12200000	pressure-resistant   pressure-resistant to 10 bar   300 l/h 950 l/h 1000 l/h   300 l/h 950 l/h 1000 l/h   1500 l 2000 l 2800 l 4000 l   0.1 – 20 µS/cm 30°C 30°C 30°C   240 x 410 240 x 490 240 x 570 240 x 700   R ¾ " R ¾ " 14 kg 18 kg 24 kg 27 kg   12115000 12200000 12280000 12400000 12400000

On request, the cartridges can also be equipped with quick-connect couplings.

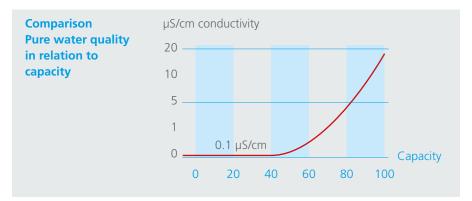

stakpure ion exchangers deliver reproducible demineralised water with a permanently low level of conductivity that remains more or less constant throughout cartridge life and meets all common standards, such as ASTM, CLSI and DIN EN 285/13060.



## stakpure industrial ion exchange cartridges

Specifically for industrial applications. With a particularly high capacity and flow rates of up to 3000 litres an hour. Made entirely of stainless steel.

stakpure mixed-bed ion exchangers	DS 7000	DS 11000	DS 15000
Water demineraliser	pressure-resistant		
Stainless-steel cartridge	pressure-resistant to 10 bar		
Flow rates	2000 l/h	2500 l/h	3000 l/h
Capacity for 10° total dissolved solids	7000 l	11000 l	15000 l
Pure water quality	0.1–20 µS/cm		
Max. water temperature	30°C		
Material	Stainless steel – AISI 316		
Dimensions mm (Ø x height)	363 x 600	363 x 850	363 x 1100
Connection	R 3⁄4 "		
Weight, approx.	55 kg	70 kg	90 kg
Order no.	12700000	12110000	12150000

On request, the cartridges can also be equipped with quick-connect couplings.

# Everything under control. stakpure digital and analogue conductivity monitoring.

### stakpure DC 400

Developed for online measurement of pure and ultrapure water downstream of ionexchange cartridges and for monitoring ring-main systems. High-precision conductivity meter ensuring exact measurement of temperature and conductivity for constantly monitoring quality. Digital display guarantees to show readings even at the lowest end of the conductivity measurement range.

Comes complete with connecting cable, plug-type power supply unit, set of two 1.5 m hoses and measuring cell.

#### Voltage selection

Automatic voltage selection for international use.

#### **LED indicators**

Show red/green to signalise when limit values are exceeded for temperature and conductivity.

Technical specifications	DC 400	
Conductivity measurement range	0.1–300 µS/cm (automatically switches over at 9.99)	
Temperature measurement range	0.1-50°C	
Temperature compensation	automatic	
Conductivity limiting-value setting	0.1–30 µS/cm, infinitely variable	
Temperature limiting-value setting	1–50°C infinitely variable	
Signals on exceeding limiting value	red/green LED indicator + buzzer (both limiting values can be deactivated)	
Fault signalling	Display changes from green to red	
Interface (optional only)	potential-free contact	
Power supply	110/230 V/50-60 Hz (automatic voltage selection)	
Dimensions mm (W x D x H)	73 x 32 x 141	
Installation	Wall mounting	
Connection	R 3⁄4 "	

Unit	DC 400	DC 400	DC 400
Version		with potential-free contact	with potential-free contact + solenoid valve
Order no.	14180500	14180600	14180700

On request, the conductivity meters can also be equipped with quick-connect couplings.

#### Large graphic display

2-colour, back-lit graphic digital display indicating conductivity in  $\mu$ S/cm or in M $\Omega$ cm and temperature in °C. Background illumination changes from green to red to signal a fault and show plaintext message.



#### Interfaces

Direct connection of peripheral equipment at potential-free contact.

#### **Limiting values**

Infinite adjustment of limiting values for temperature and conductivity.

#### **High-precision measurement**

Automatic calibration before each measurement with integrated reference resistor, temperature measured by platinum temperature sensor.

### stakpure AC 100

Conductivity meter as the easy and reliable answer to measuring the conductivity of pure water. Analogue display of measurement value in  $\mu$ S/cm. Either for connecting directly to ion-exchange cartridges or for space-saving wall installation.

Comes complete with measuring cell and set of two 1.5 m hoses.

Unit	AC 100	AW 100
Measurement range	0.1–50 µS/cm	0.1–50 µS/cm
Power supply	230 V/50-60 Hz	230 V/50-60 Hz
Dimensions mm (W x D x H)	93 x 123 x 170	93 x 125 x 96
Connection	R 3⁄4″	R ¾″
Order no. (for connecting directly to DS 1500–15000)	14160100	
Order no. (for wall installation only)		14160200

On request, the conductivity meters can also be equipped with quick-connect couplings.



AC 100



# stakpure

**stakpure GmbH** Auf dem Kesseling 11 D 56414 Niederahr Phone: +49 (0) 2602 10673-0 Fax: +49 (0) 2602 10673-200 info@stakpure.de **www.stakpure.de** 





We are certified according to ISO 9001: 2015

Is reliable and economic preparation of pure and/or ultrapure water a topic for you? Just call us!

info@stakpure.de www.stakpure.de

Retailer panel

Subject to change and printing errors.